

MSG 175A - RED FD12 BDS-05 PROCEDURE UPDATE

1
2
3
4
5
6
7
8
9
10
11
12
13
14
15
16
17
18
19
20
21
22
23
24
25
26
27
28
29
30
31
32
33
34
35
36
37
38
39
40
41
42
43
44
45
46
47
48
49
50
51

Once BDS-05 BZK Wipes have been depleted (approx. MET 11/06:00), Alcohol Wipes can be used in their place for the remainder of the mission. For your reference, this affects the procedures in 5 places on the MEDIA SAMPLE AND PCBA ANALYSIS CUE CARD (BDS-05 Experiment Ops Checklist):

1. In Step 1, delete 'BZK Wipes (5)'
2. Add the following new step after step 2:
Caddy 2 2a. Unstow Alcohol Wipes (3) from Alcohol Wipe Kit, temp stow Wipes
3. Add to Step 10 as follows:
Was: 10. Don and clean gloves
Is: 10. Don and clean gloves **using an Alcohol Wipe**
4. Add the following bold text to Step 11:
Was: 11. Swab OUTLET PORT
Is: 11. Swab OUTLET PORT **using an Alcohol Wipe**
5. Change Step 38 as follows:
Was: 38. Swab OUTLET PORT and Media Sample Bag Septum with BZK
Is: 38. Swab OUTLET PORT and Media Sample Bag Septum with **Alcohol Wipe**